


Instadose®VUE Wireless Dosimeter

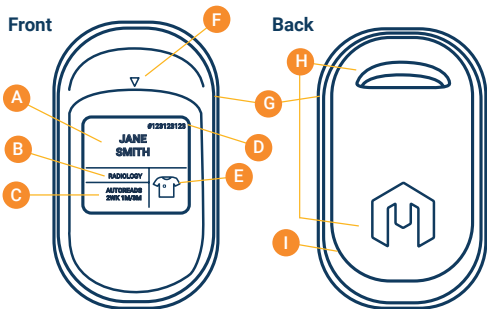

- | | |
|---|--|
| A Wearer's Name | F Detector Location |
| B Location / Department | G On-Demand Read Button |
| C Auto-Read Schedule | H Clip/Lanyard Holder |
| D Account Number | I Serial Number
(Located Under Clip) |
| E Badge Wear Location
(Body Region) | |

Dosimeter Status and Feedback Icons



Collar Badge



Torso Badge



Fetal Badge



Area Badge



Establishing
Connection
(On-Demand Reads)



Uploading
Dose Data
(On-Demand Reads)



Successful
Communication
(On-Demand Reads)



47+ Days Since
Last Successful
Communication



Unsuccessful
Communication
(On-Demand Reads)



Unsuccessful
Communication
(Calendar-set Reads)



Call
Support Team



Recall
Initiated



Badge Error:
Low Temp.



Badge Error:
High Temp.



Badge Error:
Fatal Temp.



Mirion Technologies, Inc.

PRODUCT NAME: Instadose
PRODUCT MODEL: VUE
FCC ID: 2AAZN-INSTAVUE
IC: 21847-INSTAVUE

This device complies with part 15 of the FCC Rules.
Operation is subject to the following two conditions:
(1) This device may not cause harmful interference, and
(2) this device must accept any interference received,
including interference that may cause undesired operation.



Instadose®VUE Wireless Dosimeter Wear, Care, and Communication Device Ranges



Do wear badge uncovered & facing out



Do not allow dosimeter to overheat.



Do not allow dosimeter to freeze.



Do not wear badge inside or covered



Do not immerse dosimeter in water.



Do not store in a car, outdoors, or in a place with extreme temperatures.



InstaLink 3 Gateway
30-foot range



Instadose Companion
5-foot range

Instadose Companion Mobile App Quick Start Steps

1. Download free app from Apple or Google Play Store.
2. Login with username & password email to you from dsd-support@mirion.com
3. Activate Bluetooth® setting on your smart device.



Scan to Download
Instadose Companion
Mobile App

Additional resources available at:
www.instadose.com/instadoseVUE

Performing On-Demand Dose Readings *

1



Press and hold Read Button for no more than 2 seconds.

2



The hourglass icon indicates device is searching for a connection.

3



The upload cloud icon indicates dose data transmission.

4



The checkmark icon indicates successful transmission.

* Automatic readings need no action.

The dosimeter will read automatically at the scheduled time if it's near a communication device.

NOTE Icons with exclamation point indicate unsuccessful transmission. Retry on-demand read for successful data transmission.



NOTE To confirm dose transmission, review dose history on your read history screen on the mobile app or online within the Account Management Portal (AMP+).